

SEIKA DIGITAL IMAGE

Date:10/18/2019

No: Sample : Deodorant Spray

Sample Information

Nozzle Name:[Average]

Style:STD

Category:

Delivery Co.:

Delivery Date:2019-10-18

Remark:n:1.36

Testing Information

Instrument:311XP

Test Range:0.1-100

Environment:

Medium:

Measuring Man:

Measuring Time:2019-10-18

Analyse Result (Analyse-Mode:Independent)

Xv10=2.510 μm

Xv50=10.766 μm

Xv90=21.490 μm

X[4,3]=11.488 μm

Xn10=0.476 μm

Xn50=1.211 μm

Xn90=2.457 μm

Mass Con : 17.5g/m³


S/V=9758.106 cm²/cm³

Number Con: 147425 /cm³

Fit Error:0.006

Optics Concentration:10.82462ncentration:0.001750%

(Note: V stands for cumulative volume distribution, N stands for Cumulative Quantity Distribution)

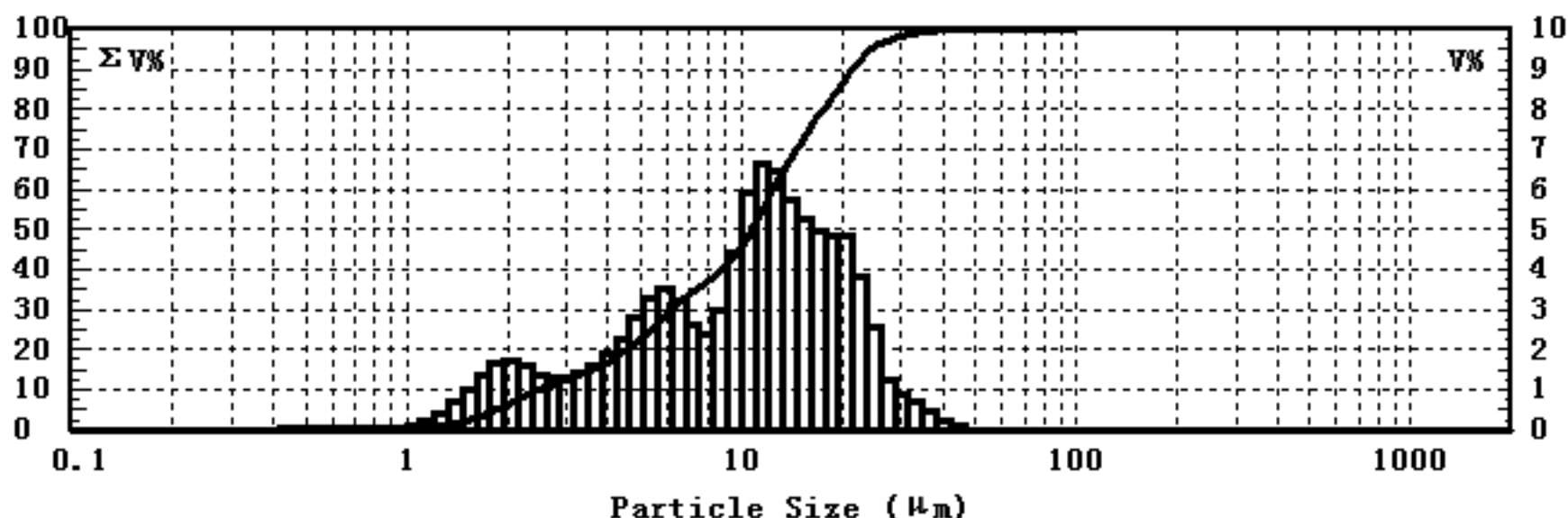

d (μm)	V%	ΣV%	d (μm)	V%	ΣV%	d (μm)	V%	ΣV%
0.383	0.000	0.000	1.957	1.670	5.893	10.000	4.404	45.476
0.422	0.008	0.008	2.154	1.731	7.624	11.007	5.947	51.423
0.464	0.018	0.026	2.371	1.581	9.205	12.115	6.622	58.045
0.511	0.033	0.059	2.610	1.365	10.570	13.335	6.458	64.503
0.562	0.048	0.107	2.873	1.271	11.842	14.678	5.769	70.271
0.619	0.057	0.164	3.162	1.292	13.134	16.156	5.278	75.549
0.681	0.057	0.221	3.481	1.399	14.533	17.783	4.935	80.484
0.750	0.051	0.272	3.831	1.594	16.127	19.573	4.833	85.318
0.825	0.045	0.318	4.217	1.884	18.011	21.544	4.814	90.132
0.909	0.046	0.364	4.642	2.286	20.296	23.714	3.807	93.939
1.000	0.064	0.427	5.109	2.775	23.071	26.102	2.553	96.492
1.101	0.113	0.540	5.623	3.274	26.345	28.730	1.218	97.710
1.212	0.223	0.764	6.190	3.509	29.854	31.623	0.892	98.601
1.334	0.412	1.176	6.813	3.237	33.091	34.807	0.671	99.272
1.468	0.678	1.853	7.499	2.636	35.727	38.312	0.430	99.702
1.616	1.001	2.855	8.254	2.354	38.081	42.170	0.221	99.923
1.778	1.368	4.223	9.085	2.991	41.072	46.416	0.077	100.000

Contact us

Name:Seika Digital Image

Tel:03-3405-1280

Address:info@seika-di.com

URL:www.seika-di.com

Printing Time:10/18/2019